

# OPERATION MANUAL

## 1. Specification

Ultrasonic Cleaner

Model : KSU- 100 / 200

Capacity : 1.6 Liter / 140x135x100(H)mm  
3.3 Liter / 240x140x100(H)mm

Frequency : BLT 40KHz

Material : Bath – Stainless Steel  
Case – ABS

Power source : AC220V, 50/60Hz

## 2. Operation

- 2.1 Fulfill the water or detergent up to 2/3 of bath.
- 2.2 Turn on the switch.
- 2.3 Turn on the timer
- 2.4 Input the materials to clean into the bath.
- 2.5 Take out the material by tools.
- 2.6 Turn off the switch after working.
- 2.7. Empty the bath in case the water / detergent were contaminated.

## 3. Attention

- 3.1 Turn on the switch in only case the water is in the bath.
- 3.2 Fulfill the cleanser the temp. is below 60oC.
- 3.3 Do not use the acid cleanser.
- 3.4 Set it at the place of well ventilated and less humidity.
- 3.5 Handle no shock against ultrasonic cleaner.

## 4. Remedy

- 4.1 Check the fuse.
- 4.2 Check the disconnection to heater cable.

“Thanks for choosing our product.”

HumanLab Instrument Co. ([www.humanlab.co.kr](http://www.humanlab.co.kr))

Office : B-401, Jaeun Bldg, #417-33, Younghwa-dong, Jangan-gu, Suwon 440821, Korea  
Phone : (82-31-2563403) / Fax : (82-31-2563404) / Email : [sales@humanlab.co.kr](mailto:sales@humanlab.co.kr)